

## Finally, true, trouble-free quick-connect hardware for HPLC!

What's new from the innovators at Optimize Technologies? Opti-Lynx™! The world's only low dispersion, high pressure quick connect system for HPLC is now available from S.I.S.

Now you can make and break High Pressure connections in seconds or reconfigure your HPLC system without tools. A quarter turn is all you need to assure leak-free connections to your analytical column, in-line filters, guard column or anywhere you frequently change components.

### Here's how it works:

**1** Start by connecting your existing tubing to an OPTI-LYNX Holder Cap using the included fingertight fitting.



**2** Then, tighten a slotted Holder Tube into the Holder cap. There are two Holder Tubes of different lengths to choose from—one for OPTI-LYNX guard column cartridges, and one for filtration elements and straight-through (union) connections.



**3** For every plug, there's a socket, and for every OPTI-LYNX Holder Tube, there's a Bayonet Connector. We have Bayonets for making ZDV connections to any analytical column (our Universal Port Adapter) and Bayonets that adapt to 10-32 fingertights, for use in-line anywhere on your HPLC system.




**4** When it's time to bring the two halves of OPTI-LYNX together for a high-pressure, quick-connect seal, all you need is an OPTI-LYNX insert. Choose a guard column cartridge, a precolumn or in-line filter, or a sealing element for straight-through connection.




## OPTI-LYNX Column Connector Kit

When a guard column is not required, OPTI-LYNX Column Connector Kit brings quick-connect convenience to any 10-32 analytical column. Use one Column Connector Kit at the top of the column for fast and easy connection with integral pre-column filtration, and another at the bottom of the column with a sealing element for a straight-through connection to the detector cell.

The basic holder set is the same whether or not you require filtration. Simply add the correct two letter extension to the part number below — CD or CE for PEEK-encased stainless steel filters, or AB for Sealing Elements.

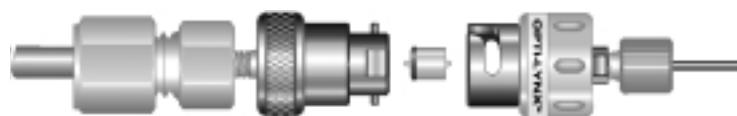

Opti-Lynx™ Column Connector Kit  
with Sealing Element

Part No.	Description	Price
11-02226AB	 Opti-Lynx Connector Kit Includes: Universal Port Adapter, Holder Cap/Holder Tube, 1 Sealing Element	